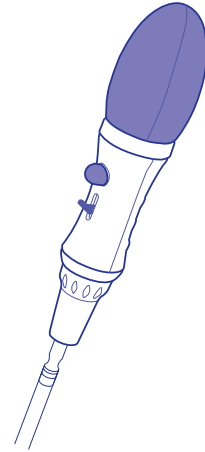
PROFILLER™ 437

MANUAL PIPETTE CONTROLLER MANUELLE PIPETTIERHILFE PIPETEUR MANUEL

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SAFETY PRECAUTIONS

- Read operating instructions carefully.
- Observe manufacturer's instructions.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals to clean instrument.
- Do not use Profiller™ 437 pipette controllers with liquids incompatible with silicone, PP, PTFE and ABS copolymer.
- Do not allow liquids to enter the instruments.

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DESCRIPTION

Profiler™ 437 is an easy-to-use manual pipette controller, which facilitates aspiration and distribution of a variety of reagents. It is compatible with glass or plastic pipettes with nominal volumes ranging from 1 to 100mL and perfectly suitable for left or right handed operation.

OPERATION

Pipette fitting: Carefully attach the pipette by inserting it into the autoclavable silicone holder of the nose cone without using force.

Aspiration: Hold instrument in a vertical position with the pipette tip down. Squeeze suction bulb. Place tip of the pipette into the liquid to be aspirated. Press lever control button slowly upward allowing the liquid to fill the pipette until the meniscus level is slightly above the required mark. The suction bulb can be squeezed again if additional aspiration is needed during pipetting. Adjust meniscus position by pressing down the lever control button.

Dispensing: Place pipette against the wall of the receiving vessel and press down lever control button to dispense.

"To deliver" pipette: when meniscus has reached a still stand, wait for the required period of time indicated on the pipette (i.e. EX +10s) for dispensing to be completed.

"Blow out" pipette: as soon as the meniscus has reached a still stand, press discharge bulb to expel the rest of the liquid contained in the pipette.

Remove the pipette: Hold the pipette and rotate gently pulling the pipette from the silicone pipette holder.

TIGHTNESS

Take care of the instrument to keep it in perfect working condition. To test tightness, attach a pipette and fill it to half of its nominal capacity. Drops falling from the pipette tip indicate a leak. Unscrew lower assembly and check silicone pipette holder and membrane filter. Reassemble correctly.

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MAINTENANCE

Do not allow liquid to enter the Profiller™ 437. The PTFE membrane of the protection filter would clog and suction be impaired. To replace the filter, remove the pipette, unscrew nose cone, pull out and replace membrane filter. The imprinted face should be toward the pipette adapter. Reassemble the unit. Clean the instrument with a damp cloth. Profiller™ 437 is not autoclavable, except silicone pipette holder.

Note: Only use the manufacturer's recommended replacement filter.

WARRANTY

The Profiller™ 437 is guaranteed for a one-year period against any factory or material defects. Damages due to non-respect of manufacturer's instructions are excluded from the warranty. Return the instrument to the dealer from whom it was purchased, after obtaining return authorisation, should regular maintenance not eliminate a detected defect.

Decontaminate instrument prior to returning it.

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